

gibco



**Cellvival
of the fittest**

Increase neuronal cell survival by more than 50% with the new Gibco B-27 Plus Neuronal Culture System

ThermoFisher
SCIENTIFIC

The standard in neuronal cell culture just got better

Since 1993, Gibco™ B-27™ Supplement and Neurobasal™ Medium has been the trusted standard for a variety of neuronal culture applications, with citations in more than 11,000 publications. However, as the desire for more reliable and biologically relevant models has increased, so too has the necessity for a media system that can maintain and mature optimal densities of functional neurons over longer periods of time in vitro. To meet this need, we have developed the new Gibco™ B-27™ Plus Neuronal Culture System. It features an optimized formulation, upgraded manufacturing process, and more stringent quality control for raw materials and final product. These improvements enable:

- Increased neuronal survival by more than 50%—the highest in the industry
- Accelerated neurite outgrowth
- Improved electrophysiological activity and maturation of neurons

The B-27 Plus Neuronal Culture System includes:

- B-27™ Plus Supplement (50X), 10 mL
- Neurobasal™ Plus Medium, 500 mL

Seamlessly replace other neuronal cell culture systems, such as classic Neurobasal Medium with B-27 Supplement in maintenance, maturation, and differentiation protocols with no change to your current workflow.



Designed to increase neuronal survival by more than 50% compared to the classic products





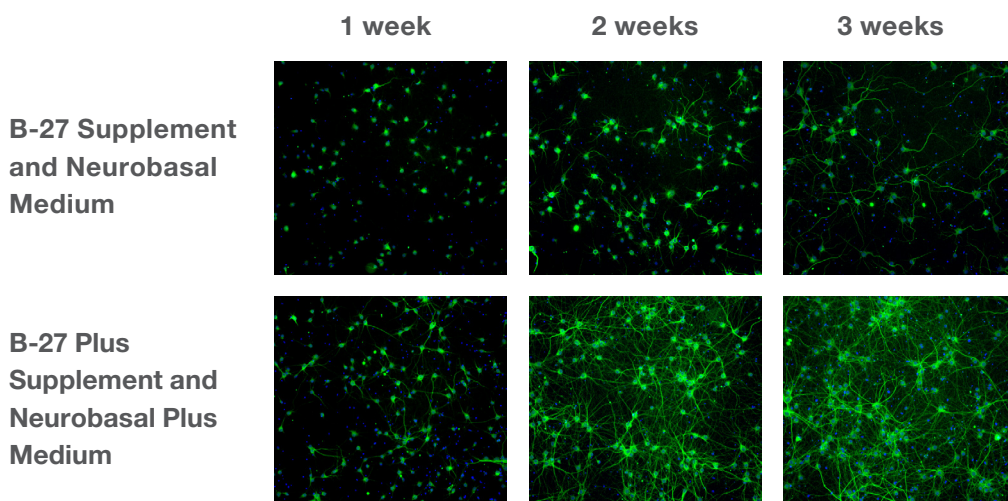






Figure 1. B-27 Plus Neuronal Culture System maintains higher neuronal survival rates compared to B-27 Supplement and Neurobasal Medium. Gibco™ Primary Rat Cortex Neurons (Cat. No. A10840) were cultured for up to 3 weeks in the listed media system and then immunostained for MAP2 on days 7, 14, and 21.

Superior culture system for primary rodent and human iPSC-derived neurons

B-27 Supplement and Neurobasal Medium

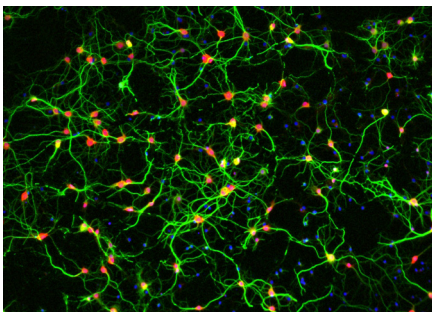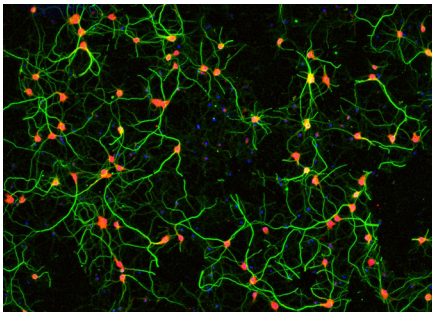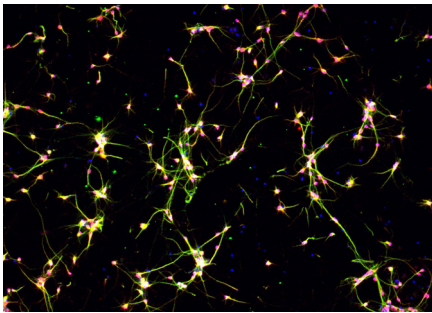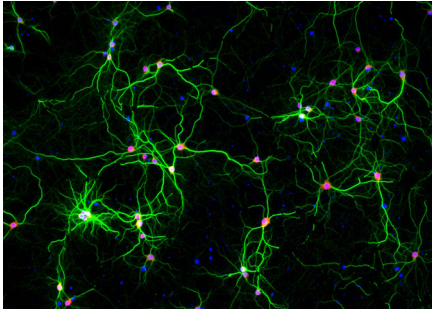

B-27 Plus Supplement and Neurobasal Plus Medium

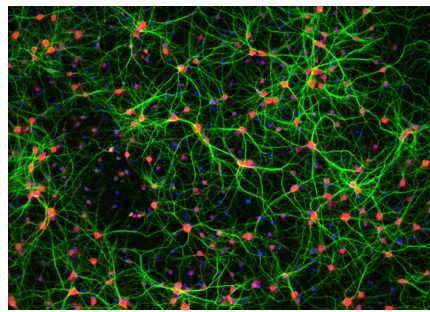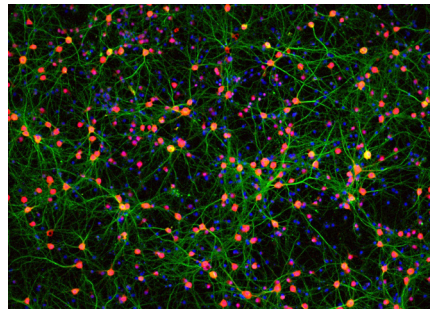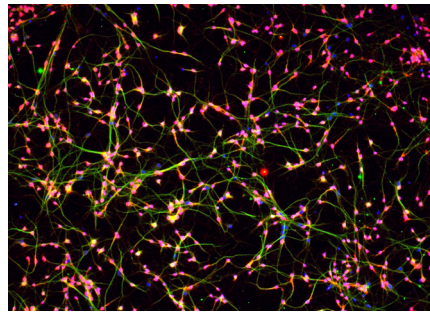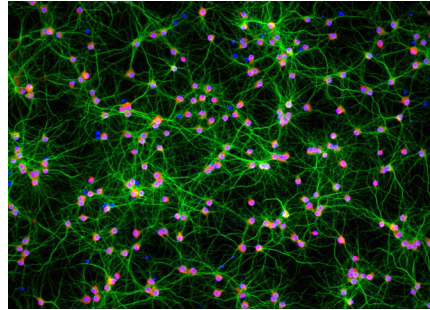

■ MAP2 ■ HuC/D ■ DAPI

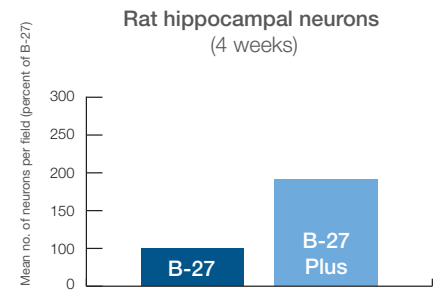
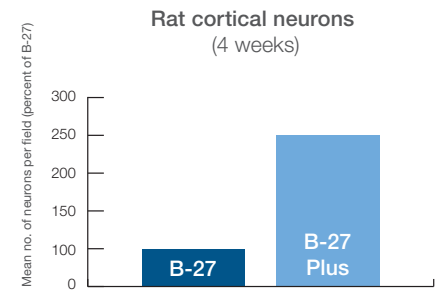
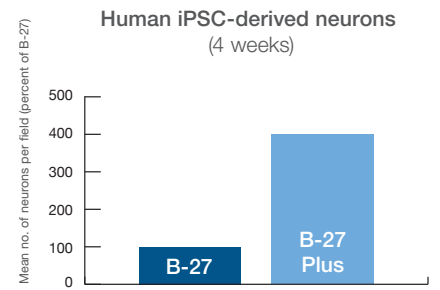
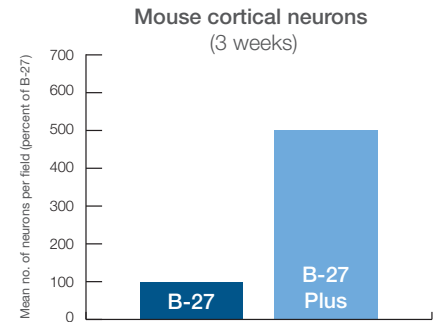


Figure 2. B-27 Plus Neuronal Culture System enables superior cultures of primary rodent and induced pluripotent stem cell (iPSC)-derived neurons compared to B-27 Supplement and Neurobasal Medium. Cryopreserved neurons were cultured for 3–4 weeks in the listed media system. Neurons were immunostained at the indicated time points for the neuronal dendritic marker MAP2 (green) and the neuronal cell body marker HuC/D (red). Nuclei were counterstained with DAPI (blue).

Supports highest survival of neurons compared to alternative media systems

Neuronal survival at three and four weeks

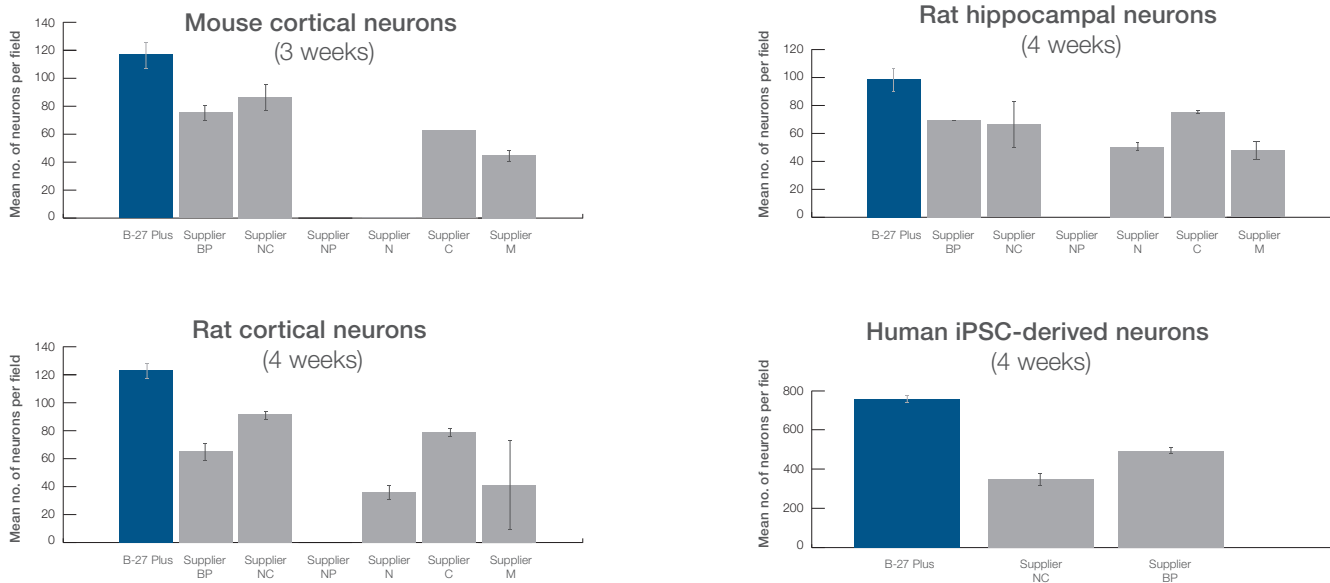


Figure 3. The B-27 Plus Neuronal Culture System achieves the highest neuronal survival in long-term cultures compared to other commercially available serum-free neuronal media systems. Primary Rat Cortex Neurons (Cat. No. A10840), Gibco™ Primary Rat Hippocampus Neurons (Cat. No. A10841), Gibco™ Primary Mouse Cortical Neurons (Cat. No. A15586), and human iPSC-derived neurons that had been matured for 9 days before cryopreservation (HIP™ Neurons, MTI-GlobalStem), were thawed in classic Neurobasal Medium with B-27 Supplement. The neurons were plated onto poly-D-lysine-coated 96-well plates and maintained for 3–4 weeks in the B-27 Plus Neuronal Culture System and alternative serum-free supplemented media systems following the suppliers’ recommended protocols. Neuronal survival was quantitated by immunofluorescent labeling using Invitrogen™ HuC/HuD Monoclonal Antibody (Cat. No. A21271) at 3 or 4 weeks. Data shown are from one of three experiments, with each run showing that B-27 Plus Neuronal Culture System enables the highest neuronal cell survival among the alternative systems.

Accelerates and increases neurite outgrowth

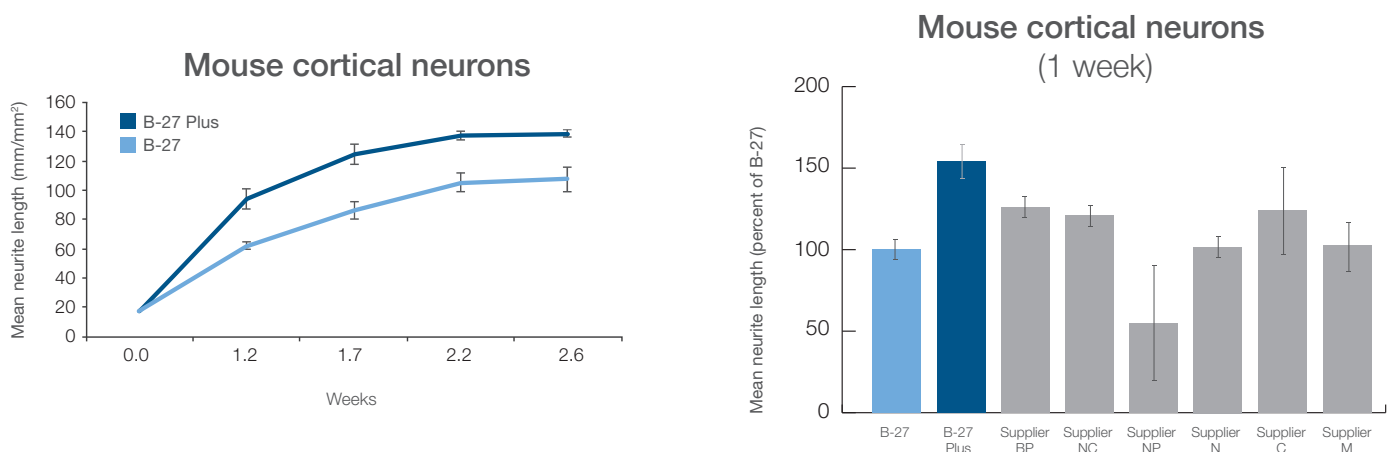


Figure 4. Acceleration of neurite outgrowth compared to B-27 Supplement and Neurobasal Medium and alternative media systems. Primary Mouse Cortical Neurons (Cat. No. A15586) were thawed in classic Neurobasal Medium with B-27 Supplement and plated onto poly-D-lysine-coated 96-well plates. Neurons were maintained for ~3 weeks in Neurobasal Medium with B-27 Supplement, Neurobasal Plus Medium with B-27 Plus Supplement, and other commercially available serum-free neuronal media systems following the suppliers’ recommended protocols. Neurite outgrowth was quantitated using differential interference contrast images taken at the time points specified. The B-27 Plus Neuronal Culture System significantly accelerates neurite outgrowth over the first few weeks compared to classic Neurobasal Medium with B-27 Supplement. Data shown are from one of three experiments, with each run showing that after 1 week, the B-27 Plus Neuronal Culture System enabled the highest mean neurite length compared to all alternative systems.

Improves electrophysiological activity and maturity


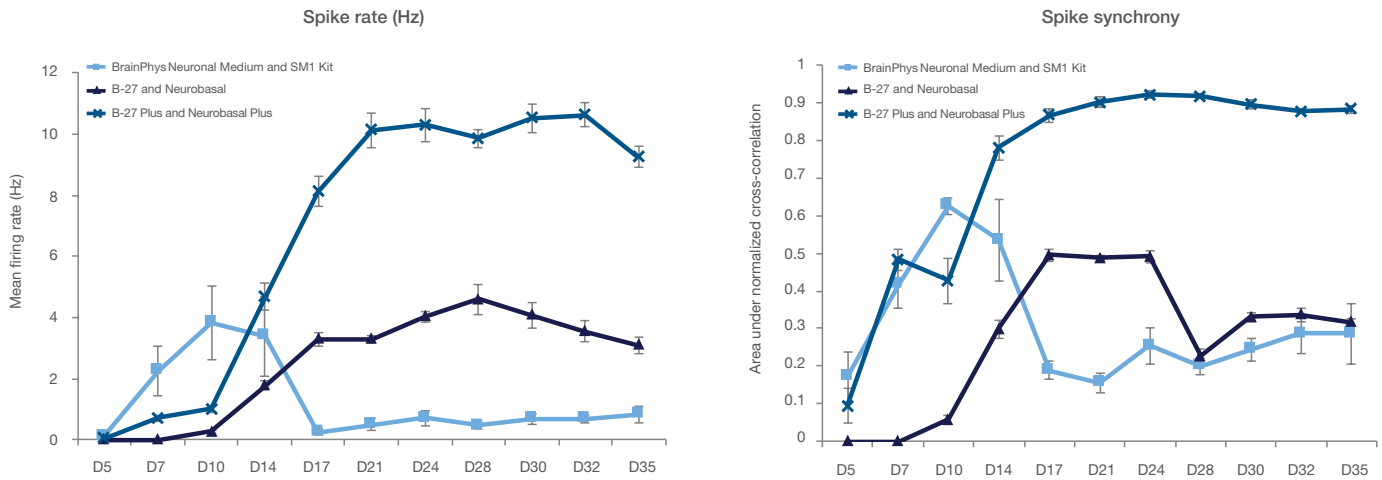


Figure 5. Improvement of electrophysiological activity compared to BrainPhys Neuronal Medium and SM1 Kit. Primary Rat Cortex Neurons (Cat. No. A10840) were plated on to 48-well multi-electrode array (MEA) plates. Cells were cultured for 35 days in the listed media system following the suppliers' recommended protocols. Spontaneous electrophysiological activity was recorded throughout with the Axion BioSystems™ Maestro™ MEA platform. Data shown are from one of four experiments, with each run showing similar trends. The B-27 Plus Neuronal Culture System enabled more consistent, stable, and highly synchronized spontaneous electrophysiological activity over time.

B-27 Supplement and Neurobasal Medium

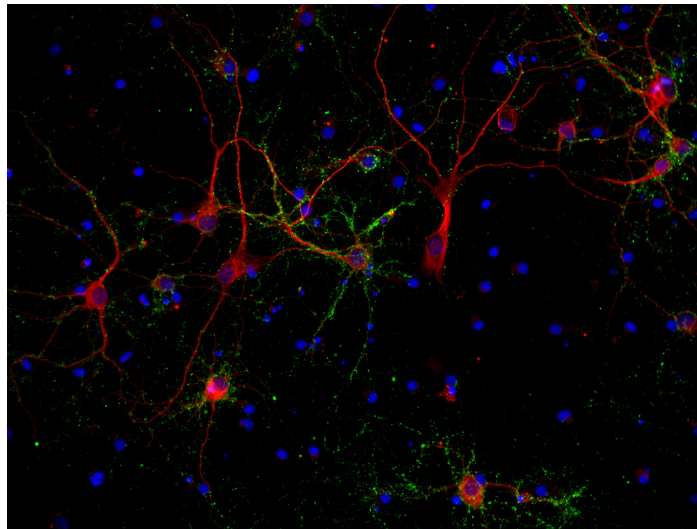

B-27 Plus Supplement and Neurobasal Plus Medium

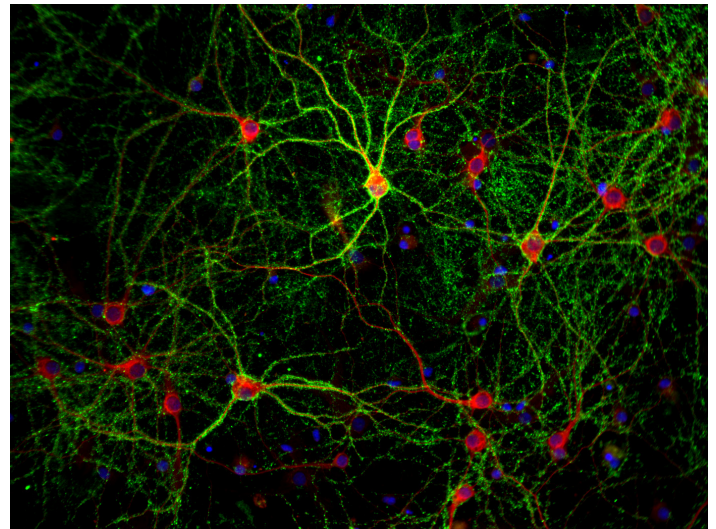


Figure 6. Enhancement of neuronal maturation compared to B-27 Supplement and Neurobasal Medium. Primary Rat Cortex Neurons (Cat. No. A10840) at day 22 were stained for the dendritic marker MAP2 (red) and synapsin 1/2 to label presynaptic terminals (green). Neurons maintained in the B-27 Plus Neuronal Culture System had significantly higher synaptic-positive puncta.

Recommended applications and ordering information

B-27 Plus Supplement and Neurobasal Plus Medium come in the same formats and will work in exactly the same protocols as the current catalog products. The classic B-27 and Neurobasal products will continue to

be offered; however, we recommend using the Plus versions and other combinations of medium and supplement in the following applications:

Application	Recommended supplement(s)	Recommended basal medium
Maintenance and maturation of prenatal primary neurons	B-27 Plus Supplement	Neurobasal Plus Medium
Differentiation, maintenance, and maturation of stem cell-derived neurons	B-27 Plus Supplement and CultureOne Supplement	Neurobasal Plus Medium
Electrophysiology studies	B-27 Plus Supplement	Neurobasal Plus Medium
Expansion of neural stem cells	StemPro NSC SFM Supplement	KnockOut DMEM/F-12 Basal Medium
Maintenance/maturation of postnatal and adult brain neurons	B-27 Supplement	Neurobasal-A Medium
Studies of insulin secretion or insulin receptors	B-27 Supplement Minus Insulin	Neurobasal-A Medium
Translational cell therapy	CTS B-27 Supplement, XenoFree*	CTS Neurobasal Medium* or CTS Neurobasal-Medium
Studies of oxidative stress and damage, apoptosis, or where free-radical damage to neurons occurs	B-27 Supplement minus antioxidants	Neurobasal Medium

Product	Quantity	Cat. No.
B-27 Plus Neuronal Culture System		A3653401
B-27 Plus Supplement (50X)	10 mL	A3582801
Neurobasal Plus Medium	500 mL	A3582901



Find out more at thermofisher.com/b27plus

ThermoFisher
SCIENTIFIC

For Research Use Only. Not for use in diagnostic procedures. © 2017 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. B-27 is a trademark of the Board of Trustees of Southern Illinois University. BrainPhys is a trademark of the Salk Institute for Biological Studies. **COL04934 0917**

* For Research Use or Manufacturing of Cell, Gene, or Tissue-Based Products. CAUTION: Not intended for direct administration into humans or animals.