detectors





Thermo Scientific[™] Dionex[™] products are UHPLC compatible by design, establishing the new standard in conventional LC. Integrating hardware, software, and separation chemistry, These products offer UHPLC to everyone for all needs.

The Thermo Scientific Dionex FLD-3000 is a high-sensitivity Fluorescence Detector series for Thermo Scientific Dionex *UltiMate*[™] 3000 HPLC systems. It is available in Rapid Separation (RS) and Standard (SD) versions. The optics of the Dionex FLD-3000 series provide maximum stray-light suppression for unrivaled detection sensitivity. Operated with the Thermo Scientific Dionex Chromeleon[™] Chromatography Data System (CDS) software, the detector provides automated qualification, various tools for method development, and instrument wellness monitoring for ease of use, maximum uptime, and the highest degree of regulatory compliance.

- Data collection at up to 200 Hz for optimal support of even the fastest UHPLC separations (Dionex FLD-3400RS).
- Standard detectors which operate at up to 100 Hz data rate for optimum support of 62 MPa (9,000 psi) Dionex UltiMate 3000 standard systems.
- Lowest limits of detection with a Raman signal-to-noise ratio (S/N):
 > 550 ASTM (> 2100 using dark signal as noise reference).
- Unsurpassed reproducibility with active flow cell temperature control for stable fluorophore activity independent of changes in ambient temperature.
- Long-life xenon flash lamp for highest sensitivity and long-term operation without the need for frequent lamp changing.



Part of Thermo Fisher Scientific

- Optional second photomultiplier (PMT) for unique Dual-PMT operation, offering an extended wavelength range up to 900 nm without sacrificing sensitivity in the standard wavelength range.
- Two-dimensional (2D) or threedimensional (3D) excitation, emission or synchro scans to provide the highest degree of flexibility for method development or routine sample characterization.
- Innovative Variable Emission Filter for real-time compoundrelated sensitivity optimization (Dionex FLD-3400RS only).
- Large front-panel display for easy monitoring of the detector status.
- Two flow-cell sizes for easy optimization to application requirements (the 8 μ L flow cell is ideal for trace analysis, and the 2 μ L flow cell offers best peak resolution with narrow-bore HPLC and UHPLC columns).

Dionex FLD-3000 Series Features	Dionex FLD-3000 Series Benefits
Innovative optical design for maximum suppression of stray light	Allows unmatched sensitivity for lowest detection limits.
Unique fast-switching Variable Emission Filter for stray light reduction in front of the emission monochromator	Generates increased detection sensitivity by adapting the filter setting to the selected emission wavelength or by recording all light emitted with wavelengths higher than the selected cutoff filter in a single data channel (Dionex FLD-3400RS systems only).
Unique Dual-PMT option	Combines flexibility with sensitivity and maximizes the wavelength range up to 900 nm without sacrificing sensitivity in the UV-vis spectral range.
Data collection rate of up to 200 Hz in single-channel mode (Dionex FLD-3400RS under Chromeleon 7.1 software control)	Provides full support of any UHPLC application. Sufficient data points even for peak widths in the subsecond range, delivering superior quantitation accuracy in ultrafast separations.
Temperature-controlled flow cells	Supplies unsurpassed reproducibility in run-to-run and day-to-day routine operation, even at fluctuating column or ambient temperature.
High-intensity xenon flash lamp with smart control features	Provides significantly longer lamp life (> 15,000 h in long life mode) than conventional, permanently illuminating xenon lamps. In addition, the lamp only consumes 10% of the power of a conventional lamp.
2D and 3D scans of emission and excitation wavelengths	Identifies quickly the optimum emission and excitation wavelength settings for your application.
8 μ L analytical flow cell 2 μ L micro flow cell	The 8 μ L analytical flow cell is the best choice for trace detection in combination with 2 mm to 4.6 mm i.d. UHPLC columns. The 2 μ L micro flow cell provides the best chromatographic efficiency and resolution with 2 mm i.d. columns. All relevant flow cell details are tracked by the detector and added to each sample.
Flow cell syringe injection kit	Cost-effective upgrade to introduce sample into the flow cell and perform an off-line measurement. Easy option to flush the flow cell when preparing for idle times.
Multiwavelength operation (Dionex FLD-3400RS systems only) on up to four channels with independent excitation and emission wavelengths	Allows detection of analytes with different fluorophores in a single analytical run.
Effective lamp age counter normalizes the lamp operation time and alerts users when to change the lamp. Various internal checks continuously monitor correct operation of all relevant detector components.	Allows the user to plan ahead for required maintenance to ensure correct operation.
Internal wavelength calibration procedures; qualification period is saved in the module	Avoids the use of unqualified equipment, even when the module is switched to another system.

Unmatched Performance Through Innovative Optical Design

The Dionex FLD-3000 series combines advanced optics with several unique features for extraordinary low-light scatter and superior S/N performance (Figure 1). The Variable Emission Filter ensures the highest possible sensitivity over the whole wavelength range. High-precision grating monochromators support fastest excitation and emission wavelength changes suitable also for multichannel operation (Dionex FLD-3400RS only). The Dual-PMT option increases the detection range up to 900 nm without the usual loss in UV sensitivity. The temperature-controlled flow cell increases the detection precision even under fluctuating ambient temperature.

A New Dimension of Sensitivity

In fluorescence detection, undesired stray light significantly increases noise, and at the same time can give a false reading for the signal value. This can make calculation of the true S/N value difficult. The optics of the Dionex FLD-3000 series are highly optimized with regards to stray light suppression. This reduces noise, ensures a true reading for the signal value, and maximizes the S/N value across the entire wavelength range (Figure 2).

Optimized Fluidics for Best UHPLC Compatibility

Dionex FLD-3000 flow cells are designed for best UHPLC support. Both the micro and analytical flow cell variants are optimized for minimum dispersion and best peak symmetry. Figure 3 shows a separation of five polycyclic aromatic hyrdrocarbons (PAHs) on a Thermo Scientific Acclaim[™] RSLC 2.1 mm × 50 mm, $3 \,\mu \text{m}$ column, detected with a $2 \,\mu \text{L}$ (micro) and an 8 μ L (analytical) flow cell. Although all peaks elute within a time window of only 0.7 min and peak volumes are only 27 μ L to 46 μ L, even the analytical flow cell detects symmetric peaks and fully resolves

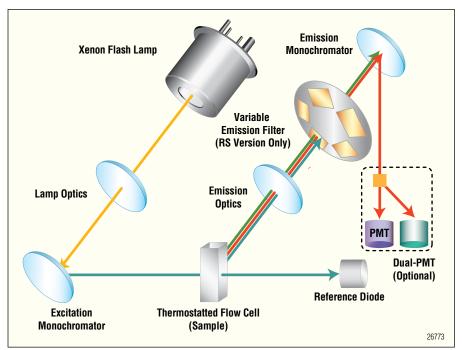


Figure 1. Schematic of the optical design of the Dionex FLD-3000 series.

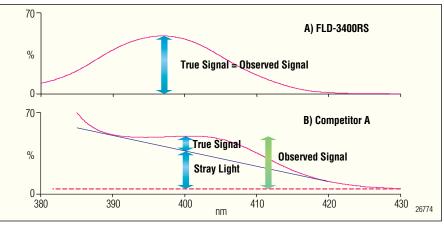


Figure 2. Raman band of water (397 nm): A) fully resolved from stray light with the Dionex FLD-3400RS; and B) poorly resolved with a competitior's fluorescence detector.

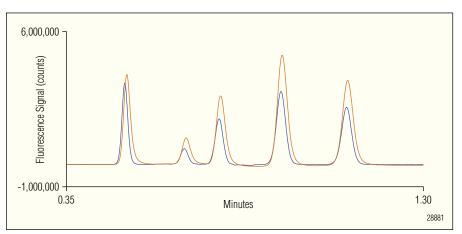


Figure 3. Sub 1.2 min separation of five PAHs on a 2.1 \times 50 mm, 3 μ m column using an analytical (orange) and a micro (blue) flow cell. The analytical flow cell resolves all peaks and achieves the best S/N. The micro flow cell provides the best resolution.

them. The best S/N is obtained with the analytical flow cell. The micro flow cell achieves the best chromatographic resolution.

RSLC Technology for Any Application

The Dionex FLD-3400RS is the technology leader in fluorescence detection and ideally matches the performance and flexibility of the RSLC platform. Data rates up to 200 Hz, very high sensitivity, tunable emission filters, two flow-cell variants, different lamp operation modes, and detection at up to 900 nm without compromise provide more possibilities than any other fluorescence detector on the market. Run existing applications or easily develop new ultrasensitive, ultrafast applications with sub-2 min run time (Figure 4).

Shortest Run Times Even with Excitation and Emission Switching

For the best fluorescence sensitivity, trace compounds must be analyzed with optimum excitation and emission wavelengths. In a separation of different fluorescing analytes, it becomes necessary to switch the wavelengths between the peaks. UHPLC conditions shorten the available time window for the required grating movements significantly. The Dionex FLD-3000 series is designed with fast-turning grating monochromators which reach the new setting in less than 250 ms. Figure 5 indicates the total switching time of 350 ms under real-life conditions, including response time and software operation delay. This time includes a sensitivity change and a response time delay. This switching time is compared against a third-party detector operated under identical conditions. The competitive instrument requires 6.1 s to reach the new settings and is therefore not suitable to support UHPLC separations. In contrast, the Dionex FLD-3000 detectors are ideally suited to support very complex UHPLC separations. Figure 6 shows a 5.5 min separation of 17 fluorescing PAHs (with eight changes of wavelength, sensitivity, and filter).

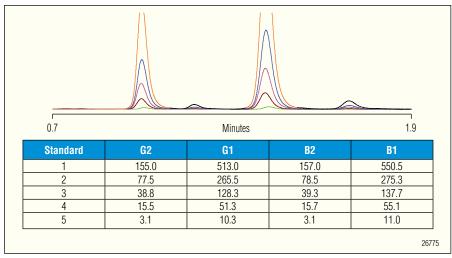


Figure 4. Overlay of five aflatoxin standard separations using an Acclaim RSLC C18 2.1 × 100 mm, 2.2 µm column (B2, B1, G2, G1 in order of elution). Table concentrations in ppb.

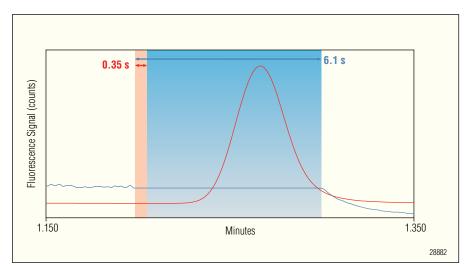


Figure 5. Data acquisition with a Dionex FLD-3000 series detector (red) and a third-party detector (blue). The Dionex FLD-3000 series detector easily completes the wavelength and sensitivity change before the peak. The third-party detector is ready after the peak has already left the flow cell.

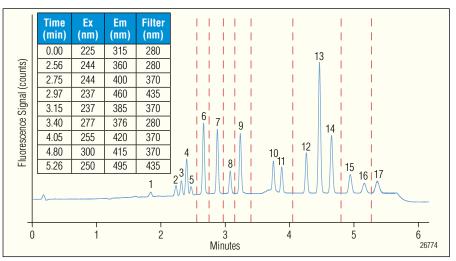


Figure 6. Fluorescence detection of 17 PAHs within 5 min using a Nucleodur C18 PAH 3×100 mm, 3 µm column at a flow of 2 mL/min. The detection is made possible by the fast-moving filters and monochromators of the Dionex FLD-3400RS.

Unsurpassed Reproducibility

Fluorescence is a function of temperature. Changing temperatures will yield different emission intensities for equal fluorophore concentrations. To minimize this effect, the flow cells of the Dionex FLD-3000 series are temperature controlled. This significantly improves the precision in run-to-run and day-to-day routine operation, even at fluctuating ambient temperatures (Figure 7).

Features for Maximum Uptime

Dionex FLD-3000 detectors use a xenon flash lamp. This lamp features three different flash frequency modes (LongLife, Standard, HighPower) for optimizing the sensitivity or lamp lifetime. Operating the lamp in LongLife mode leads to lifetimes of >15,000 h, several times more than a conventional, continuouslyilluminating xenon lamp used in other manufacturers' detectors. Conventional lamps also require a significant warmup period and must be turned on even if no fluorescence signal is acquired. Therefore, the real data acquisition time of a conventional xenon lamp can be significantly lower than its specified lifetime. The Dionex FLD-3000 series flash lamp is thermally optimized for warm-up within seconds. This means the lamp is used only when it is needed-during the acquisition.

Smart Use of Lamp Frequency Settings

The different lamp frequencies can be varied even during a chromatographic separation. Combining the lamp frequencies in a smart way provides optimum S/N and increases lamp lifetime at the same time. During dwell times and equilibration, the lamp flash frequency can be reduced (LongLife, 20 Hz). While peaks are eluting, a highlamp-flash frequency (e.g., HighPower, 300 Hz) reduces baseline noise and, therefore, results in optimum S/N. In the Figure 9 example, the reduction of the flash frequency during idle times increases the lamp lifetime by 100%.

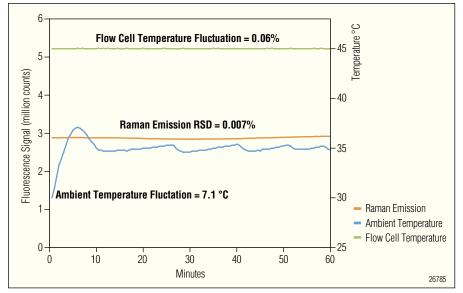


Figure 7. The thermostatted flow cell guarantees reproducible results even under varying ambient temperatures. Overlay of Raman signal intensity (orange), flow cell temperature (green) and ambient temperature (blue).

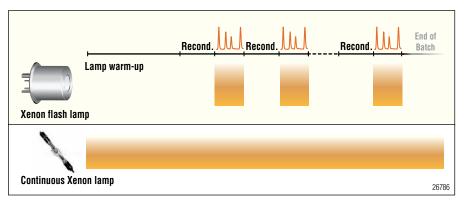


Figure 8. Lamp handling for maximum lifetime. The xenon flash lamp does not require long warm-up times and is only on during data acquisition. For minimal lamp wear, it can even be switched off when no peaks elute, e.g., during reconditioning times. Continuous xenon lamps are less flexible and require long-term temperature stabilization..

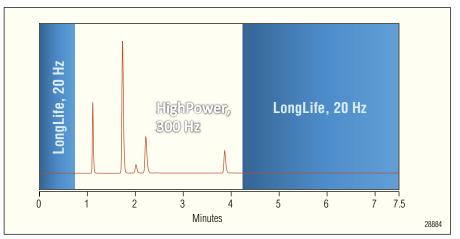


Figure 9. Alternating lamp flash frequencies during a run combines best S/N with an increased lamp lifetime. In this example, reducing the number of lamp flashes during idle times increases lamp lifetime by 100%.

Combining Sensitivity with Flexibility

PMTs are commonly used for fluorescence detection due to their significantly higher sensitivity compared to photodiodes. However, PMTs provide a limited wavelength range. A typical PMT supports an emission wavelength range up to 600 nm. To overcome this limitation, the Dionex FLD-3000 detectors can be ordered with, or field upgraded to, a Dual-PMT configuration. This unique option adds a second red-sensitive PMT to extend the wavelength range to 900 nm without sacrificing the sensitivity of the UV-vis region (Figure 10). Even the PMT selection can be changed during a chromatographic run to make sure each analyte is detected at highest S/N. Figure 11 provides an example of a dye separation. In addition to the switching of the excitation and the emission wavelengths, the detecting PMT changes whenever beneficial.

The Ultimate in Flexible Method Development

The Dionex FLD-3000 series provide various tools for effective development of selective methods for fluorescent analytes.

- 2D or 3D Excitation, Emission, or Synchro Scans: Perform a scan at a defined time of the chromatogram or simply use the 3D scanning functionality. This mode repeatedly acquires spectra during the entire separation and allows reviewing the data with familiar DAD-like spectral tools (Figure 12).
- Unique Variable Emission Filter: This automatically picks the best suitable emission filter for maximum stray light reduction (Dionex FLD-3400RS only).
- Zero-Order Mode: When the fluorescence characteristics of an analyte are unknown, the zero-order mode of the monochromator can be used to optimize the separation. It sets the emission grating to reflect the complete emission spectrum on the PMT.
- Multichannel Operation: Simultaneous operation of up to four different excitation and emission wavelengths are possible (Dionex FLD-3400RS only).

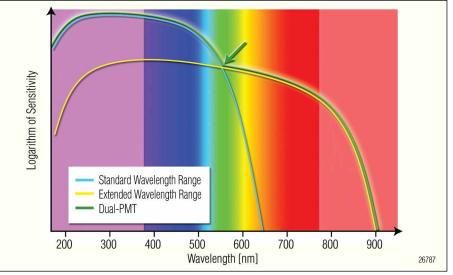


Figure 10. The Dual-PMT option of the Dionex FLD-3000 series is a unique solution for providing optimum sensitivity across all wavelengths.

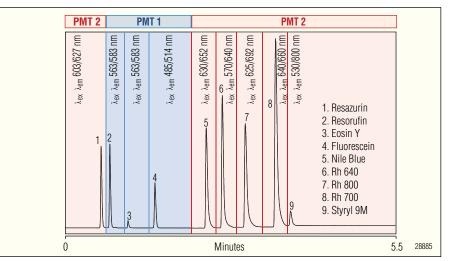


Figure 11. UHPLC analysis of dyes with a Dual-PMT detector configuration, which can change the measuring PMT during a chromatographic run. This ensures that any analyte achieves optimum S/N across the entire wavelength range up to 900 nm.

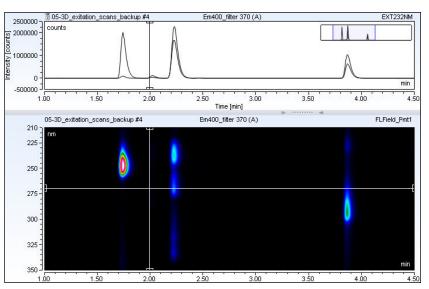


Figure 12. Excitation, emission, and synchro 3D scans acquire spectral data during the entire chromatographic run. From a single run, optimum detection wavelengths can be extracted by interfacing with an easy-to-use, DAD-like 3D field console.

Take Control of All Functions—with Ease

Using Operational Simplicity[™] as its guiding principle, the latest generation of Chromeleon software, version 7.1, takes users from samples to results in the shortest time possible. This software suite ensures users are utilizing the Dionex FLD-3000 detectors effectively and with ease, despite its various options and functions. In addition to the intuitive instrument control shown in Figure 13, Chromeleon's main new processing features include the Cobra[™] peak detection algorithm for fast, accurate peak detection, and the SmartPeaks[™] Integration Assistant for easy handling of unresolved peaks. Users can also automate processes with eWorkflows, a new tool that automates all chromatography workflows via an intuitive and easy-to-use interface.

Thermo Scientific Dionex AutoQ Equipment Qualification

Dionex AutoQ[™] Equipment Qualification routines turn tedious qualification work into a simple task. The Chromeleon CDS automatically operates the following test protocols:

- Installation Qualification
- Operational Qualification
- Performance Qualification

Upon completion, the software creates comprehensive reports, including passed/failed checks and charts.

Method Transfer Made Easy

Conventional LC methods can be made much more productive by using short UHPLC columns packed with smaller particles. The Dionex UltiMate 3000 systems fully support any UHPLC column format; the only remaining challenge is to define the right application parameters. Chromeleon 7.1 helps in the process of transferring a method to a different column format. It features an embedded Method Transfer Wizard that combines chromatographic results already available in the CDS, with some user input on the planned column to create a new instrument method (Figure 15).

Based on the prediction of the chromatographic performance, the user can decide to further speed up the method or calculate for a different column. The Wizard also predicts how much time and solvent the new method will save.

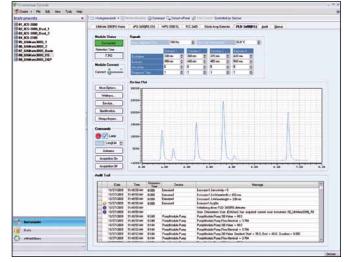


Figure 13. Intuitive ePanels provide direct access to all main instrument control features and easy access to advanced options.

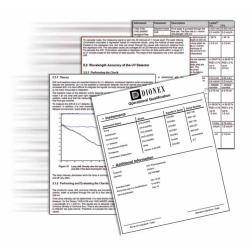


Figure 14. Operational qualification documentation for the Dionex UltiMate 3000 System.

(9) (0. (d. /* here	SHELO	luthod, Transfer (Leap	ince) - Ownedian Ohi	mangaphy Dude			
and the states	Reference of Houses	je kont finge i e	(e bury Constant) (e bury Constant) (e bury Constant)	27	and I and East and East	U	
Optiment Method *	Vers Joges Conversion AdvancedCore Conversion Parameters				canced Conversel	Method Executor	_
O Derrore	Please specify the driver To further accelerate se			e new odkew.			U
6 (PE SHOWS)		Current Colum	n Now Column				
Frencher fullP1-3000	Longth	250.0 9	100.0	(10.0.1000.0 and			
L Gardon	Dismeter	45	30 💌	\$1.1 . 100 0 mm]			
PAG-MORES	Paricle sce	50 .	25 .	(0.1			
See para-success) See fasteri	O Boost factor		1.18	+ 0.68 million			
(h lino	· Flow	0.000	8.773	(0.0018.000 wil/in	ie)		
	Pressure limit		1034	101034 bad			
HPLC_Instance(_method HPLC_restrict_2 CHPLC_restrict_2	* Results	Current Colu	nn New Column	s Saving	9		
injection List	Resolution factor (EP)		1.31				
instrument Method	Max perceae Inaction volume	200 ber 12.000 ut	727 bar 2 283 µl		#1 X		
/v Data Processing	Elverturage	30.750 mi	5.232 ml		83.5		
2 Report Designer	Runtime	28 500 mm	1510 min	1510 min 22 %			
Electronic Report	Thoughput		+5.7				
Spectral Library							

Figure 15.Transfer of conventional LC methods is exceptionally easy with the Method Transfer Wizard embedded into Chromeleon 7.1.

KEY SPECIFICATIONS

Light Source:				Excitation/Emission Wavelength Switching Time:		
Xenon Flash Lamp				<250 ms		
Lamp Pulse Frequency:				Wavelength Accuracy:		
HighPower (~300 Hz); Standard (~100 Hz); LongLife (~20 Hz)			ngLife (~	±2 nm		
Data Collection Rate:				Wavelength Repeatability:		
Single wavelength: up to 200 Hz (FLD-3400RS under Chromeleon 7.1 control)			under	±0.2 nm		
Multiple wavelength: up to 4 Hz (FLD-3400RS))	Sensitivity:		
Single wavelength: up to 100 Hz (FLD-3100)				Raman S/N: > 550 ASTM over the entire lifetime of the		
Spectrum Scanning Modes:					lamp (> 2100 using dark signal as noise reference)	
2D or 3D excitation, emission or synchro scans (3D scans under			(3D scan	Flow Cell Thermostatting:		
Chromeleon 7.1 control)				Ambient + 15 °C to 50 °C		
Wavelength Range:					Flow Cells:	
Variant	Ex Min	Ex Max	Em Min	Em Max	Analytical flow cell volume: 8 μ L	
FLD-3100	200	630	265	650	20 bar (290 psi) maximum cell pressure	
FLD-3400RS	200	630	220	650	Micro flow cell volume: $2 \mu L$	
FLD-3100 with Dual-PMT	200	880	265	900	40 bar (580 psi) maximum cell pressure	
FLD-3400RS with Dual-PM7	200	880	220	900	PC Connection:	
Spectral Bandwidth:					All functions controllable through USB	
Excitation: 20 nm					GLP Features:	
Emission: 20 nm					Automatic Equipment Qualification (Dionex AutoQ), System	
Emission Filter:					Wellness, and Qualification Monitoring with Chromeleon,	
Variable, 5 positions (FLD-34	00RS)				cell i.d. chip	
Fixed, 280 nm (FLD-3100)					Dimensions $(h \times w \times d)$:	
No. of Fluorescence Channels:					$16 \times 42 \times 51 \text{ cm} (6.3 \times 16.5 \times 20 \text{ in.})$	
Up to 4 (FLD-3400RS)					Weight:	
1 (FLD-3100)					18 kg	

ORDERING INFORMATION

In the U.S. and Canada, call 1-847-295-7500, or contact the Dionex Regional Office nearest you. Outside of North America, order through your local Dionex office or distributor. Refer to the following part numbers.

Detectors

Dionex FLD-3100 Fluorescence Detector (without flow cell)	
Dionex FLD-3100 Fluorescence Detector with Dual-PMT (without flow cell)	
Dionex FLD-3400RS Rapid Separation Fluorescence Detector (without flow cell)	
Dionex FLD-3400RS Rapid Separation Fluorescence Detector with Dual-PMT (without flow cell)	
Accessories	
Analytical Flow Cell, 8 μ L, stainless steel	
Micro Flow Cell, 2 µL, stainless steel	
Dual-PMT Option	
Dual 1 M1 Option	
Flow Cell Syringe Injection Kit	

Nucleodur is a registered trademark of MACHEREY-NAGEL GMBH & CO. All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsideries.

Nucleodur is a registered trademark of MACHERET-NAGEL GMBH & CO. All other trademarks are the property of Thermo Pisner Scientific Inc. and its subsider

Speed · Simplicity · Solutions

North America

Dionex Products

1228 Titan Way P.O. Box 3603 Sunnyvale, CA 94088-3603 (408) 737-0700

E

U.S./Canada (847) 295-7500 South America Brazil (55) 11 3731 5140

Europe

Austria (43) 1 616 51 25 Benelux (31) 20 683 9768; (32) 3 353 4294 Denmark (45) 36 36 90 90 France (33) 1 39 30 01 10 Germany (49) 6126 991 0 Ireland (353) 1 644 0064 Italy (39) 02 51 62 1267 Sweden (46) 8 473 3380 Switzerland (41) 62 205 9966 United Kingdom (44) 1276 691722

Asia Pacific

Australia (61) 2 9420 5233 China (852) 2428 3282 India (91) 22 2764 2735 Japan (81) 6 6885 1213 Korea (82) 2 2653 2580 Singapore (65) 6289 1190 Taiwan (886) 2 8751 66551



www.thermoscientific.com/dionex

LPN 2381-02 PDF 0811 ©2011 Thermo Fisher Scientific, Inc.

DIONEX