


EASY-nLC 1200 System

Best-in-class nano LC, preconfigured for straightforward state-of-the-art separations

The Thermo Scientific™ EASY-nLC™ 1200 system provides effortless peak performance for nanoflow applications up to 1200 bar. Dual in-line flow sensors before solvent mixing give excellent gradient precision. Its intuitive software facilitates rapid method setup and immediate system control.

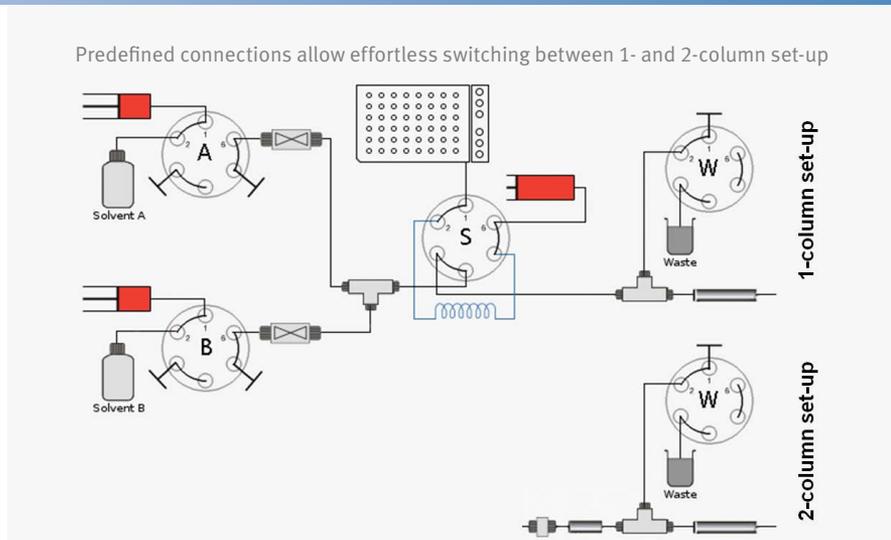

Increased System Robustness and Servicability

- New maintenance-free ceramic valves result in extraordinary dependability and lower cost of ownership
- Thermo Scientific™ nanoViper™ connections for tool-free, reliable and fast connections
- Easy access to all serviceable parts

Intelligent Maintenance

- Automated, built-in maintenance routines
- User-friendly leak tests pinpoint position of leaks
- Advanced pressure test quickly diagnoses blocked capillaries
- Detailed event log for troubleshooting
- Remote access capability for system interrogation and technical support

Improved Performance and Throughput at up to 1200 bar

- Faster sample loading and column equilibration increase sample throughput without sacrificing performance
- Use of longer analytical columns for higher peak capacity

Advanced Flow Control


- Intelligent Flow Control (IFC™) during loading and equilibration steps provides robust performance while increasing throughput
- Automatic Flow Control (AFC™) during gradient delivery provides outstanding reproducibility throughout injection sequence



Product Specification


Integrated EASY-nLC Software and Touch Screen

Direct system control of the EASY-nLC system via touch screen interface allows for fast system set up, method configuration, and maintenance.

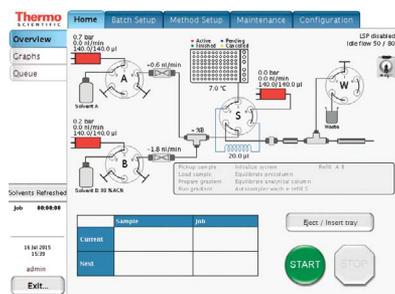

Xcalibur Data System

Thermo Scientific™ Xcalibur™ software is the state-of-the-art solution for controlling Thermo Scientific mass spectrometers. The EASY-nLC software is fully integrated with Xcalibur software, allowing full LC-MS control through a single software solution.

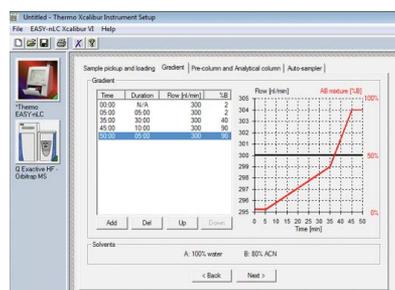

MS Interfaces

The EASY-nLC 1200 system was designed to work together with Thermo Scientific™ EASY-Spray™ Ion Sources and Nanospray Flex™ Ion Sources, making it compatible with EASY-Spray columns and all other nano LC columns.

Easy MS Connection

EASY-nLC instruments integrate seamlessly with Thermo Scientific mass spectrometers, including:

- Thermo Scientific™ Q Exactive™ series instruments
- Thermo Scientific™ LTQ Orbitrap™ series instruments
- Thermo Scientific™ Orbitrap Fusion™ Tribrid™ instruments
- Thermo Scientific™ TSQ™ Triple Quad instruments

For more information, visit: thermofisher.com/nanoLCMS

©2016 Thermo Fisher Scientific Inc. All rights reserved. PEEK is a trademark of Vitrex PLC. All other trademarks are the property of Thermo Fisher Scientific and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Africa +43 1 333 50 34 0
Australia +61 3 9757 4300
Austria +43 810 282 206
Belgium +32 53 73 42 41
Canada +1 800 530 8447
China 800 810 5118 (free call domestic)
 400 650 5118

Denmark +45 70 23 62 60
Europe-Other +43 1 333 50 34 0
Finland +358 10 3292 200
France +33 1 60 92 48 00
Germany +49 6103 408 1014
India +91 22 6742 9494
Italy +39 02 950 591

Japan +81 45 453 9100
Korea +82 2 3420 8600
Latin America +1 561 688 8700
Middle East +43 1 333 50 34 0
Netherlands +31 76 579 55 55
New Zealand +64 9 980 6700
Norway +46 8 556 468 00

Russia/CIS +43 1 333 50 34 0
Singapore +65 6289 1190
Spain +34 914 845 965
Sweden +46 8 556 468 00
Switzerland +41 61 716 77 00
UK +44 1442 233555
USA +1 800 532 4752

PRODUCT SPECIFICATIONS

Specification	Value
Pressure Range	0–1200 bar
Flow Range (Gradient)	20–2000 nL/min Recommended: 100–1000 nL/min
Flow Range (Loading & Equilibration)	≤25 µL/min
Retention Time Reproducibility	0.1–0.4% RSD within recommended flow range
Pick-up Volume Range	0.10–18.00 µL (20 µL loop)
Injection Reproducibility	≤0.2% RSD at 5 µL pick-up ≤3.0% RSD at 0.1 µL pick-up
Injection Linearity	R ≥ 0.9985 for 0.5–10 µL injection volume R ≥ 0.9995 for 0.3–1.6 µL injection volume
Carry-over	<0.05% (caffeine) <0.1% for m/z 722.3 and <1.00% in general (100 fmol BSA digest, YIC*DNQDTISSK)
Gradient Delay Volume (Mixing tee to venting tee)	<1 µL
Injection Loops	20 µL default (5 µL and 50 µL optional)
Gradient and Injection Pumps	Single action syringe pumps (sapphire piston), 140 µL volume Gradient pumps: 100 µL/min refill speed Injection pump: 40 µL/min refill speed
Valves	4 identical 6-port, 3-position proprietary valves from Thermo Fisher Scientific
Solvent Compatibility	≤95% acetonitrile in water. Wetted parts: Fused silica, PEEK™, 316 stainless steel, zirconium dioxide
Sample Tray Formats	6 × 8 HPLC vials 1 × 96 or 1 × 384 MTP and 6 vials 2 × 48 or 4 × 96 PCR strips and 6 vials
Autosampler Cooling Range	Minimum of 5 °C Maximum of 20 °C below ambient
MS Drivers	LC Devices 3.0 for XCalibur 3.0 and later
Remote Support	Direct internet access for online diagnosis by EASY-nLC technical experts
Operating Conditions	5–30 °C ambient temperature 20–80% relative humidity
Dimensions (H x W x D)	45 cm × 36 cm × 38 cm (18" × 14" × 15")
Weight	37 kg (82 lb)
Main Power Supply	100–240 V, 50–60 Hz, 250 W

Ordering Information

Description	Part Number
EASY-nLC 1200	LC140
Nanospray Flex Ion Source (NG)	ES071 (ES072)
EASY-Spray Ion Source (NG)	ES081 (ES082)

Thermo
SCIENTIFIC

A Thermo Fisher Scientific Brand